INSTRUCTION MANUAL

LCD Digital Microscope (LDM)

Model # 44342

Introduction

Thank you for purchasing the Celestron LCD Digital Microscope (LDM) with a 2" monitor. Your microscope is a precision optical instrument, made of the highest quality materials to ensure durability and long life. It is designed to give you a lifetime of pleasure with a minimal amount of maintenance.

Before attempting to use your microscope, please read through the instructions to familiarize yourself with the functions and operations to maximize your enjoyment and usage. See the microscope photos to locate the parts discussed in this manual.

This microscope provides high powers from 24x up to 240x (up to 1920x with digital zoom). This microscope is ideally suited for examining specimen slides of yeasts and molds, cultures, plant and animal parts, fibers, bacteria, etc.

The LCD Digital Microscope does not use eyepieces that are used in traditional microscopes. You will view specimens on the LCD screen which are easy to see and you can enjoy them with others also. You can also take snapshots or short videos with the built-in digital camera.

The final section provides simple care and maintenance tips for you to follow to ensure that your microscope provides you with years of quality performance, usage, and enjoyment.

Figure 1

Standard Accessories Included with your Microscope

• Digital Camera – built-in	• Slide Holder
• LCD Monitor	• AC Adapter
• 4x Objective Lens	• AC Adapter Plugs – USA/EU/UK
• 10x Objective Lens	• 5 Prepared Slides, 10 Blank Slides, and 10 Cover Slips
• 40x Objective Lens	• SD Card Slot
• Bottom Illuminator	• USB Mini Cable

Specifications

Model # 44342	Specifications
Stage	Simple with Slide Holder Mechanism - 1.7"x1.7" (44mmx44mm)
Digital Camera	2MP CMOS; 6x magnification in lieu of an eyepiece; 1600x 1200 pixel array
LCD Monitor	2" with 8x Digital Zoom – High Definition Digital TFT display 640 x 240 pixels
Focuser	Smooth, single motion
Objectives	Achromatic 4x, 10x and 40x – all optical glass
Flash Memory	32MB – approximately 65 snapshots (JPEG files) @ 2MP resolution or a few minutes of VGA video (AVI files)
USB Cable	2.0 Mini – length 6.5' (2m)
Slide Holder	Accepts Standard Glass Slides – 1" x 3" (25mm x 76mm)
Power Knob	Triple with click stop
Illuminator	Built-in Electric – LED 6Volt and 6Watt
AC Adapter	Input – Universal 100 to 240Volt, 50/60HZ
Weight/Dimensions	28oz (794g) / 6.5" (165mm) x 6.5" (165mm) x 8.5" (215mm)

Magnification (Power) Table

Use the following table to determine the magnification of the different objective lenses in combination with your microscope using the normal mode of the digital image on the LCD screen and using the digital zoom feature.

Objective Lens	4x	10x	40x	
Digital Image – normal	24x	60x	240x	
Maximum with 8x digital zoom feature	192x	480x	1920x	

Setting Up Your Microscope

1. Carefully remove the microscope and other parts from the carton and set them on a table, desk, or other flat surface.
2. Remove the plastic cover from the microscope.

Figure 2

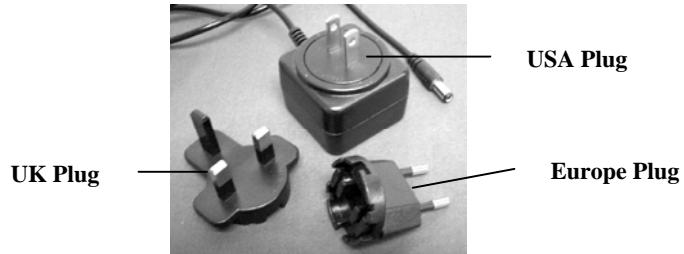
Figure 3

3. Plug the small cable from the AC adapter into the socket on the back of the base (see Figure 2 above).
4. Insert the plug end of the AC adapter into the proper power source.

Note: The AC Adapter supplied with this microscope has an interchangeable plug system that can work in the USA, UK, Europe, and many other parts of the world. The plug for the USA will be installed in the AC Adapter. You can easily change plugs for your particular requirement by doing the following:

- A. Push down on the small button labeled "PUSH" and hold it down while at the same time hold the prongs of the plug and rotate it slightly counterclockwise and pull up to remove it. Then remove your finger from holding the "PUSH" button down.
- B. Take the plug style that you want to use and center it over the AC Adapter and rotate it until it falls down into the opening. Then rotate it clockwise until you hear a click sound which lets you know the plug is now installed correctly.

Microscope Operation

Before looking at slide specimens you must remove the protective cover over the LCD screen, turn the LCD on, turn on the illuminator, and understand how to use the slide holder and then you are ready to begin viewing.

Figure 4

LCD Module

This digital microscope is different than traditional microscopes --- instead of using eyepieces to look at a specimen in a traditional microscope, the LCD monitor replaces the eyepieces so you can look at the specimen on the screen by yourself or share the views with others. To begin to view specimens with your microscope, you will have to turn the LCD monitor on by turning the Power Button “On” (see Figure 4). Push the button once and shortly the Monitor On LED (red color) will come on indicating the LCD monitor is “On”. That is basically all you need to do to use the LCD screen for viewing specimens. The various keys on the LCD Module are mainly used for taking images (snapshots and video) and will be discussed later in this manual.

Illumination

1. To turn the illuminator on, see Figure 5 below and turn the control dial to the left.
2. The illuminator is used for specimen slides where the light shines up through the hole in the stage through the slide.
3. Rotate the illuminator dial to the brightness level desired when viewing specimens.

Figure 5

Adjusting the Lighting

Specimens of different size, thickness, and color variations will require different levels of illumination. There are two ways to change the amount of illumination when viewing a specimen; adjusting the brightness on the dial shown above in Figure 5 and changing the EV function on the LCD monitor. The EV (exposure value for brightness) function increases or decreases the brightness level by using the Left and Right arrow keys on the LCD monitor.

When viewing a specimen that is not transparent or dark in color, you may need to increase the amount of light to resolve certain features or details. This is best done by simply increasing the brightness of the illuminator by rotating the brightness control dial all the way to its highest setting.

Optimum lighting will be found by experimenting with adjustments as each specimen may require slightly different illumination as well as the same specimens viewed under different powers.

Viewing a Specimen

Your instrument is provided with a fixed, simple stage with holding clips to guide the positioning of the slide holder.

1. Place a specimen slide - 1" x 3" size (25mm x 76mm) into the slide holder. Do this by gently placing the specimen slide under the small guides of the slide holder (Figure 6). Then, push the slide all the way into the cutout in the slide holder.
2. The slide holder with slide inserted will be placed onto the stage with the arrow on the slide holder pointed forward. Inside the stage (Figures 7 & 8) are two small holding clips under which the slide holder is to be inserted.
3. Push the slide holder in until you begin to see the specimen (remember to start at the lowest power). You can examine the specimen by moving the slide holder back and forth slowly and moving it slightly to the left or right.
4. For slides that have the specimen off to one side you may not be able to see the entire specimen and in these cases you can then remove the slide from the slide holder and rotate the slide 180° and try again.
5. Manually moving the slide holder may seem difficult at first but after experimenting it becomes quite easy to use.

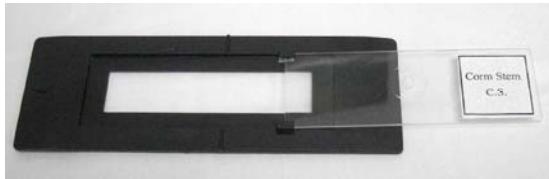
Figure 6

Figure 7

Figure 8

Once the slide is in place you will be ready to start viewing by setting the power.

6. Use the power knob (in the center of Figure 9) to rotate the objective lenses until the 4x dot is lined up near the indicator – it will click when positioned correctly. Always start with the lowest power objective (4x with this microscope) which gives you 24 power and work your way up to higher powers. At 24 power you will have the widest field of view and the brightest image.
7. Look at the LCD screen while turning the focus control (Figure 9) until the specimen comes into view. You may need to position the slide specimen as discussed above so the image is in the field of view.
8. With the 4x objective lens, you can also vary the power anywhere from 24x to 192x by using the digital zoom of the LCD module. Use the direction keys (Up/Down) on the LCD module to use the digital zoom. On the left of the screen will be an icon indicating what digital zoom setting you are at. **Note:** Digital zoom is best up to 4x.
9. For higher powers, you will need to rotate the power knob to the 10x objective position for medium power and to the 40x objective for the maximum power. You will have to refocus when changing the power of the objective lenses. While using either of the 10x or 40x lenses you also can increase power by using the digital zoom.

Figure 9

Note: At maximum powers the images may not be as sharp and dark. Your best views will be at the lower powers.

Digital Imaging

You can take snapshots or a short video with your microscope using the built-in digital camera. With the internal storage memory, you do not need to use a PC or any other devices to do imaging. Transferring of images to your PC for saving and or printing them is easy and will be discussed later in this manual.

Note: If you are going to take images, do not connect the USB cable to your PC or damages could occur to the equipment. The USB cable is not used at all for taking images!

General Settings

Before setting up the snapshot or video menus you need to make some general setups first. You can enter the general settings from the Snapshot Mode, Video Mode, or Time Mode (all three with icons showing in the top right of the LCD screen). You navigate the different modes by pressing the Esc key.

You begin by pressing the Menu key and this will have “Capture” or “Video Output” highlighted at the top left of the screen. Then press the Right direction key which will then show “Setup” highlighted. Your choices are below and you navigate to each selection by using the Up/Down keys.

1. Format – press OK to enter, choose to Cancel or Execute, and then press OK to set.
2. Language – press OK to enter, choose English, French, German, Italian, Spanish, Portuguese, simplified Chinese, traditional Chinese, or Japanese and then press OK to set.
3. Date Input – press OK to enter. Choose Off or Set. If you choose Off, then press OK to set. If choosing “Set” you will set the Day/Month/Year and time. For making the date/time settings:

- A. Use the Up/Down keys for your choice of what to display first – the day, the month, or year.
- B. Then, use the Left/Right keys to highlight the day, month, year, hour (24-hour format), minutes, or seconds. At each selection, you use the Up/Down keys to make your choices and then press OK to set the choices.
- C. Press the Left key twice to Exit the Menu.
- D. The date and time will be shown in the top right of the screen if you “Set” it.

Note: Most Settings will have to be set each time you turn on the LCD screen. You will see an “M” in the bottom right of the screen which means you are using the internal memory. If using an optional SD Card, then the SD Card icon will be shown in the bottom right of the screen.

To set up for Snapshot or Video Modes, you press the Esc key until the correct icon is in the top right of the LCD screen.

Set Up for the Snapshot Mode

The icon of a digital camera should be in the top right of the LCD screen to begin. Then press the Menu key.

1. In the top left of the LCD screen is the EV (Exposure Value) which shows the brightness level. The normal (default) position is 0.0 and this can be adjusted in increments up to +2.0 to -2.0 by pushing the Left/Right directional keys. If the default position is chosen, then the EV will not be shown on the screen.
2. In the lower right of the screen is shown the remaining snapshots available to take from the internal memory.
3. In the top right of the screen is shown an icon of a camera which indicates it is ready for snapshot mode. The resolution being used is shown in the bottom left of the screen.
4. By pushing the directional keys (Up/Down) you change the digital zoom shown by an icon.
5. Push the Menu Key to make the following choices in the highlighted “Capture” section. Push the Up/Down keys to switch to the choices and push OK to enter and set the item.
 - A. Size – this sets the resolution you want to use for the image. The camera itself uses the 1600 x 1200 pixel arrangement (2MP) sensor but you can use digital interpolation to increase the resolution or you can use lower resolutions to increase the number of images you can save in your internal storage memory. Use the directional keys to make changes and then click on “OK” to save your setting.
 - B. EV – you can change the default setting for the exposure value.
 - C. Color – you can change from Standard (normal), to Monochrome (black/white), Vivid, or Sepia.
 - D. Time Stamp – if you desire to have the time alone, date alone, or the time and date together stamped on your images you can choose which, or you can choose to have no stamp. Press OK to set your choice.
 - E. Press the Left key twice to Exit the Menu.

Set Up for the Video Mode

The icon of a video camera should be in the top right of the LCD screen to begin.

1. In the lower left of the screen is shown the resolution for taking videos.
2. In the lower right is the available time left in memory for taking videos.
3. By pushing the directional keys (Up/Down) you change the digital zoom as shown by an icon.
4. Push the Menu Key to make the following choices in the highlighted “Video Output” section. Push the Up/Down keys to switch to the choices and push OK to enter and set the item.
 - A. Size – this sets the resolution you want for the image. You should choose 800x600 pixels. Lower resolution (640x480 pixels) will allow for longer videos using the internal memory but may not be compatible with various computers. Use the Up/Down keys to make your selection and then press OK to set.
 - B. EV – you can change the default setting for the exposure value you wish to use.
 - C. Press the Left key twice to Exit the Menu.

Taking Images

1. Snapshots – to take snapshots make sure you are in the camera snapshot mode (camera icon in the top right of the screen) and you will use the Snap Key to take an image. Adjust to the sharpest focus using the focus dial as this is critical to obtaining the best images possible. Push the key when ready and the image will be stored in the internal storage memory. Experiment at best focus position and varying brightness levels and after a while you will be able to take good images.
2. Video – to take a video make sure you are in the video mode (video/movie camera icon in the top right of the screen). Use the Snap Key to begin the video and also to end it. The elapsed time of the video is shown on the screen in the bottom right.
3. You can also take images (snapshot or video) by using an SD (Secure Digital) Card (up to 1GB). Your SD card is inserted in the LCD Monitor (Figure 1). When you insert the SD card properly the SD card icon will be visible at the bottom right of the LCD screen. On the bottom right of the screen will also be shown the memory available on the SD card. When imaging this way the images are saved on the SD card and not in the internal storage memory.

Note: Inserting or removing an SD card while the LCD is on may cause the LCD to shut down and/or could damage the SD card. Always turn off the LCD screen to insert or remove an SD card.

Reviewing and Managing Your Images

You can review and manage your snapshots and videos on the LCD Monitor that are saved in the internal storage memory.

1. Push the ESC key until you enter the Stored Images Mode.
 - A. Use the directional keys to select the image you want to view – all images (and videos) will be numbered.
 - B. The selected image will have a yellow border around it and press the OK key to view this image.
 - C. When reviewing a snapshot image, you can use the digital zoom for enlargement. But, when done reviewing the image return the digital zoom to the zero position. Press the OK key to return to the stored images.
 - D. When reviewing a video image (the video/movie camera icon will be displayed on the image), you can stop and start the video by pushing the Up key. Push the Down key at any time during the video to return to the start position. For videos, in the top right of the screen is shown the length of the video and while playing the video the elapsed time is shown in the bottom center of the screen. Press the OK key to return to the stored images.
 - E. Push the ESC key to exit the stored images and return to real time.
2. When you have an image with the yellow border around it, you can manage it by pushing the Menu key:
 - A. By using the Up/Down keys you have a choice of “Delete”, “Copy To Card”, or “Protect”.
 - B. Delete – press OK to enter this selection. With the Up/Down keys you have a choice of “Single” which allows you to delete the file selected by pressing OK, “All” which allows you to delete all files in memory by pressing OK – if you select “All” you will be asked to confirm this by choosing “Execute” or “Cancel”. If you do not want to delete any files press the Left key to exit back to the Menu items.
 - C. Copy To Card – this has no function.
 - D. Protect (lock) – press OK to enter this selection. With the Up/Down keys you have a choice of “Single” which protects the image selected, “All” which protects all images in memory, or “Select” which takes you back to the stored images. For any images locked, an icon of a lock will appear at the center bottom of the screen. To “Unlock” any or all locked images, turn the LCD screen off and none of the images will be protected.
 - E. From the Menu you may push the Left key to exit the Menu.

Transferring Your Images

To transfer images to a PC, you need to have a free USB port and have an imaging program(s) for snapshots and videos.

Note: Do Not disconnect the USB cable while transferring images or damage may occur.

1. You can transfer images from the internal storage memory to your PC by using the supplied Mini USB Cable. The small plug end of the cable plugs into the LCD Monitor (Figure 1) and the large plug end of the cable plugs into your PC. Your PC will automatically recognize the new hardware. Then, you will be asked to choose which program on your PC you want to transfer the images to – you may have to choose a different program for snapshot images and another program for video images depending on what software programs are installed on your computer.
2. If you used an SD card to store your images, you can transfer them to your PC using the method in # 1 above or you can take the SD card out and use the SD card slot on your PC to transfer the images. Using either way your PC will ask you to choose which program you want to transfer the images to.




Care, Maintenance, and Warranty

Your Celestron microscope is a precision optical instrument and should be treated with care at all times. Follow these care and maintenance suggestions and your microscope will need very little maintenance throughout its lifetime.

- When you are done using your microscope, remove any specimens left on the stage.
- Turn off the illuminator switch.
- Turn off the LCD monitor.
- Unplug the power cord.
- Always place the plastic bag over the microscope when not in use or when being stored.
- Store the microscope in a dry and clean place.
- Be very careful if using your microscope in direct sun light to prevent damage to the microscope or your eyes.
- When moving your microscope, carry it with both hands (if possible) by the base or the round body.
- Clean the outside surfaces with a moist cloth.
- Always unplug any cords before cleaning.
- Never clean optical surfaces with cloth or paper towels as they can scratch optical surfaces easily.
- Blow off dust with a camel's hair brush or an air blower from optical surfaces.
- To clean fingerprints off of optical surfaces, use a lens cleaning agent and lens tissue available at most photo outlets and when cleaning do not rub in circles as this may cause sleeks or scratches to occur.
- Never disassemble or clean internal optical surfaces. This should be done by qualified technicians at the factory or other authorized repair facilities.
- When handling glass specimen slides, use care as the edges can be sharp.

Sample Images

The following specimen images left to right below were taken at 40x, 100x, 400x of a Corn Stem C.S.

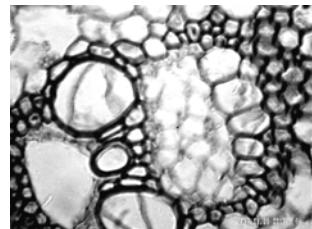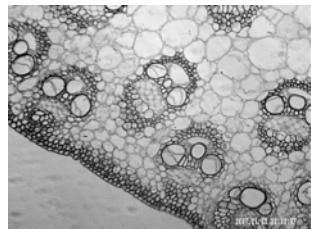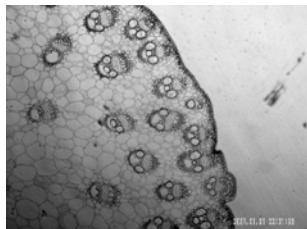Warranty

Your microscope has a two year limited warranty. Please see the Celestron website for detailed information on all Celestron microscopes at www.celestron.com.

EEC: This product complies with EEC guidelines in EN61558-2-6:1997 and EN61558-1:1997+A1

FCC Statement

This device complies with Part 15 of FCC Rules. Operation is subject to the following two conditions:

1. This device may not cause harmful interference, and
2. This device must accept any interference received, including interference that may cause undesired operation.

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Torrance, California 90503 U.S.A.